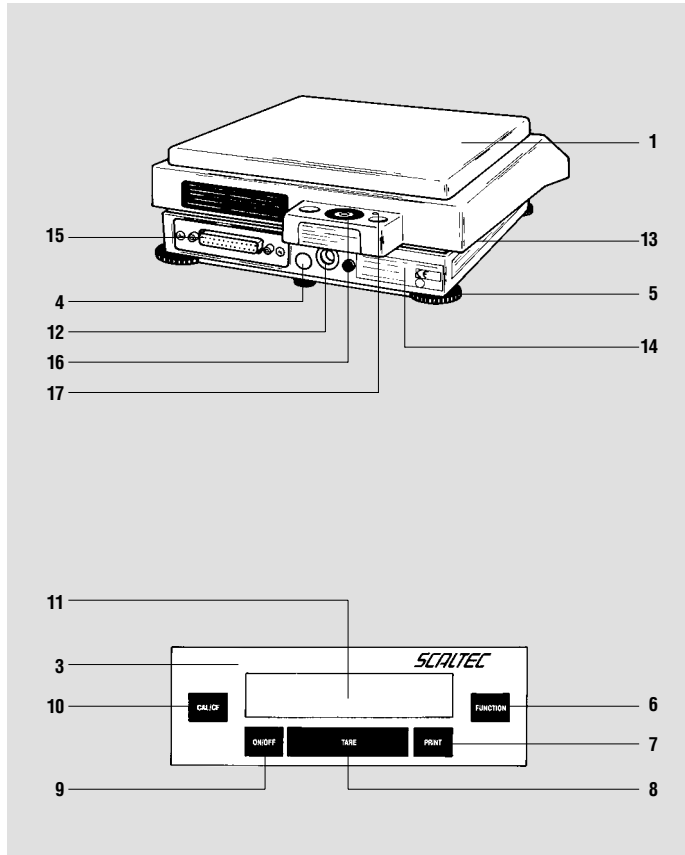
*SCALTEC*

## **SBA/SBC**

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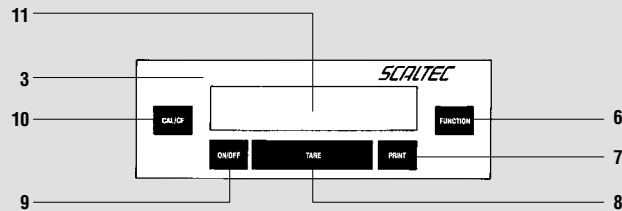
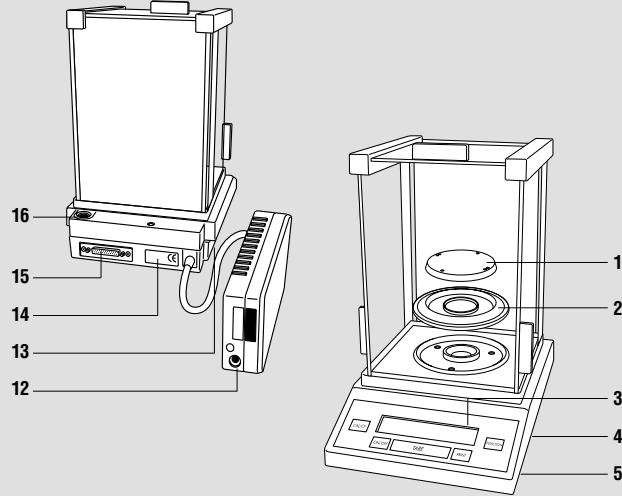
<sup>1)</sup> To activate the CF function,  
press it for more than two seconds

Pos. Bezeichnung  
No. Designation

- 10** CAL/CF-Taste<sup>1)</sup>  
CAL/CF key (calibration/clear function)<sup>1)</sup>
- 11** Gewichtsanzeige  
Weight display
- 12** Betriebsspannungsanschluss  
AC jack
- 13** Kennzeichnungsschild mit Angabe  
der eichtechnisch relevanten Daten  
Verification ID label with metrological data  
for verified balances approved  
for use as legal measuring instruments
- 14** Typenschild mit **CE**-Zeichen  
Manufacturer's ID label with the **CE** mark  
of conformity
- 15** Datenschnittstelle  
Data interface port
- 16** Libelle  
Level indicator
- 17** Befestigungsöse für Diebstahlsicherung  
Lug for fastening an antitheft locking device

# Analysenwaage Analytical Balances

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<sup>1)</sup> To activate the CF function, press it for more than two seconds

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**Before reading these instructions, fold out the first page  
showing a general view of the balance.**

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---

Please read these installation and operating instructions carefully before you begin to operate your new balance. The SCALTEC product you have purchased meets high-quality standards and is designed to assure long years of trouble-free service.

### **Warranty**

Do not miss out on the benefits of our extended warranty. Please complete the warranty registration card and return it to the address listed below (see supplement).

In the unlikely event that your balance should require servicing within the 3-year warranty period, please contact the office listed below for your country, and state the model number, serial number and the problem:

Federal Republic of Germany:  
SCALTEC INSTRUMENTS GmbH  
Robert-Bosch-Breite 10  
37079 Goettingen, Germany  
Phone (+49/551) 209773-0  
Telefax (+49/551) 209773-9  
Internet: <http://www.scaltec.com>  
E-mail: [info@scaltec.com](mailto:info@scaltec.com)

### **Storage and Shipping Conditions**

Allowable storage temperature: +5 °C ... +40 °C  
+41 °F ... +104 °F

After unpacking the balance, please check it immediately for any visible damage as a result of rough handling during shipment. If this is the case, proceed as directed in the section entitled "Safety Inspection."

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### Ambient Conditions

The SCALTEC SBA/SBC Balance is designed to provide reliable weighing results under normal ambient conditions in the laboratory and in industry.

Therefore, please observe the following when choosing a suitable place of installation:

- Set up the balance on a stable, even surface
- Protect the balance from aggressive chemical vapors
- Protect the balance from drafts that come from open windows and doors
- Avoid exposing the balance to extreme vibrations
- Avoid exposing the balance to extreme heat or to direct sunlight

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For information on the current legal requirements of your country, please contact your local SCALTEC customer service office.

### **Information on Radio Frequency Interference**

#### **Warning!**

This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause interference to radio communications. It has been tested and found to comply with the limits for a Class A computing device pursuant to Subpart J of Part 15 of FCC rules, which are designed to provide reasonable protection against such interference, when operated in a commercial environment. Operation of this equipment in a residential area is likely to cause interference, in which case the user, at his own expense, will be required to take whatever measures may be required to correct the interference.

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**Using Verified Balances  
as Legal Measuring Instruments  
You must calibrate the balance at the place  
of installation before using it as a legal measuring  
instrument** (see "Calibration," page 17).

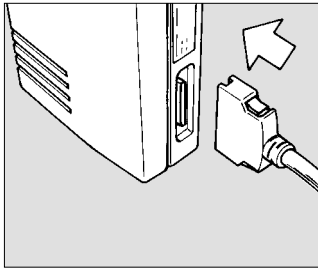
This balance is not allowed to be used for weighing goods intended for direct sale to the public. The type-approval certificate for verification applies only to non-automatic weighing instruments; for automatic operation with or without auxiliary measuring devices, you must comply with the regulations of your country applicable to the place of installation of your balance. A suitable thermometer and barometer are recommended for monitoring the ambient conditions.

For balances of accuracy class **(I)**, a thermometer and barometer are recommended for monitoring the ambient conditions. The temperature range indicated on the verification ID label must not be exceeded during operation. The balance must warm up for at least 24 hours after initial connection to AC power or after a relatively long power outage.

The manufacturer complies with EC Directive No. 90/384/EEC for non-automatic weighing instruments, which has been in effect since January 1, 1993, within the Single European Market as well as the accreditation of the quality management system of the manufacturer by Lower Saxony's Regional Administrative Department of Legal Metrology (Niedersächsische Landesverwaltungsamt-Eichwesen) from February 15, 1993.

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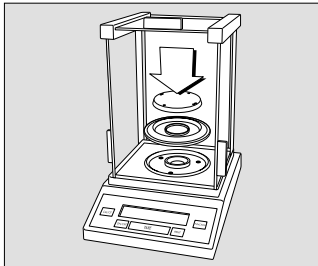
### Important Note Concerning All Verified Balances Approved for Use as Legal Measuring Instruments:

Provided that an official lead seal is required for the verified balance, a control seal is affixed to the balance. Unauthorized attempts to remove this seal will irreversibly damage it. If you break the seal, the validity of the verification will become void, and you must have your balance re-verified.

### Connecting the SBC 22 Balance to the Electronic Box:

- Plug the cable into the socket of the electronics box.

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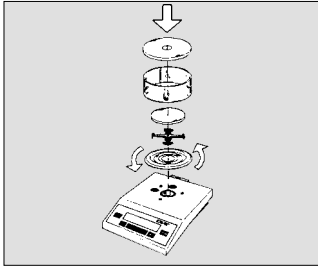
- Schirmring
- Waagschale

### Preparing the SBA 31/32/33 and the SBC 22/31/32/33 with a Draft Shield Chamber

Place the following components inside the chamber in the order given:

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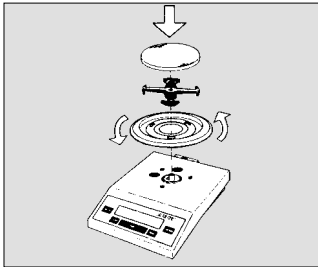
### Preparing the SBA 41/42 and the SBC 41/42 with a Round Glass Draft Shield

- Place the shield disk on the balance. Turn the disk counterclockwise until it is securely fastened.

Place components listed below on the balance in the order given:

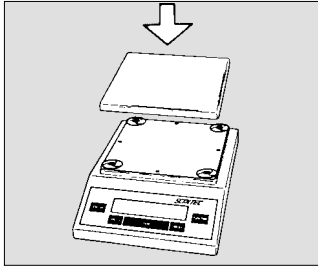
- Pan support
- Weighing pan
- Glass draft shield cylinder
- Draft shield cover

### Preparing the SBA 53 and the SBC 53 with a Round Weighing Pan

- Place the shield disk on the balance. Turn the disk counterclockwise until it is securely fastened.

Place components listed below on the balance in the order given:

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
### **Preparing the SBA 51/52/61/62 with a Rectangular Weighing Pan**

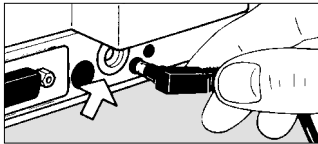– Place the weighing pan **(1)** on the balance

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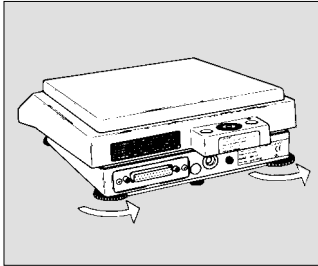
### **Connecting the Balance to AC Power**

The balance is powered by an AC adapter. Make sure that the voltage rating printed on this unit is identical to your local line voltage.

**Plug the cord of the AC adapter into the DC jack (12); then plug the AC adapter into a wall outlet (mains).**

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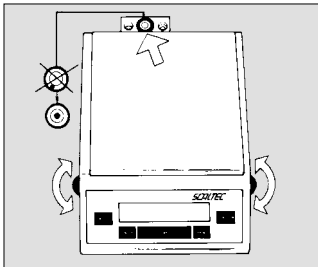
### Leveling the Balance Using the Level Indicator

At the place of installation, level the balance using the leveling feet **(5)** so that the air bubble is centered within the circle of the level indicator **(16)**.

Retract the two auxiliary feet located at the back of the balance.

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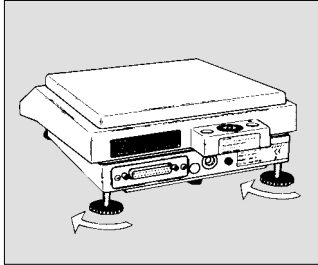
Use the level indicator as a guide to level the balance:

To lift the balance, extend the leveling feet (turn clockwise).

To lower the balance, retract the leveling feet (turn counterclockwise).

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After retracting the rear auxiliary feet, extend them until they touch the surface on which the balance rests.

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
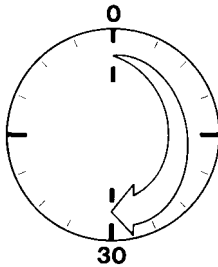
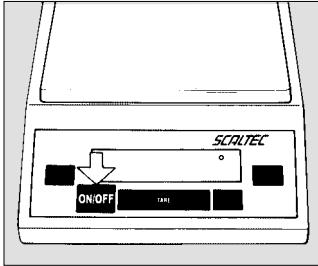
To deliver exact results, the balance must warm up for at least 30 minutes after initial connection to AC power or after a relatively long power outage. Only after this time will the balance have reached the required operating temperature.

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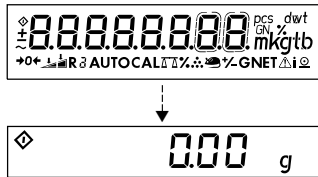### Turning the Display On or Off (Standby Mode)

Press the ON/OFF key (9) to turn the display on or off.

### Self-Test

After the balance has been turned on, an automatic self-test of the balance's electronic circuitry is performed. At the end of the self-test, a zero readout is displayed. This means that the balance is ready to operate.

### Important Note Concerning Verified Balances Approved for Use as Legal Measuring Instruments:

For verified balances that have a verification scale interval "e" which is greater than the scale interval "d," the last digit on the display is bordered.

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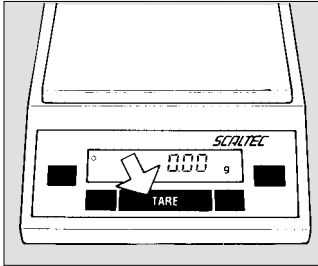
◇ für **Busy**

The display shows the following special codes for your information:

○ displayed in the upper right corner stands for **OFF**

○ displayed in the lower left corner means **standby**

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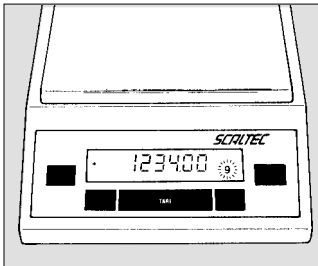
### Taring

A weight can be determined accurately only from a defined zero point. Press TARE **(8)** to zero the weight display. You can tare within the entire weighing range of the balance.

### Important Note Concerning Verified Balances Approved for Use as Legal Measuring Instruments:

The small circle in the weight display (on the left) shows that the balance is exactly tared to "0" ( $\pm 0.25$  of a scale interval).

### Simple Weighing

Place your sample on the weighing pan **(1)** to determine the weight. Read off the weight indicated on the display only after the weight unit "g" or a different unit selected appears as the stability symbol.

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During calibration, the sensitivity of the balance is adjusted to changes in the ambient conditions.

Calibrate your new balance at the place of installation after each warmup period and before the first weight measurement. In verified models the balance should be calibrated at least once a day.

**The balance offers you various calibration functions. You can select these functions by setting the appropriate menu code.**

**All SBC models have a built-in calibration weight.**

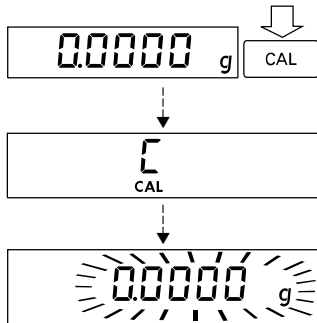
**Using Verified SBC Balances as Legal Measuring Instruments**

Before using your balance as a legal measuring instrument, you must carry out an "Internal Calibration" operation at the place of installation after the warmup period.

Menu code selection: 1 9 3 \*

The calibration weight is internally applied by servomotor and removed at the end of calibration.

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When the display shows a zero readout, press the CAL key **(10)** to activate calibration.

If any interference affects calibration, you may obtain a brief display of the error message "Err 02." In this case, tare and press the CAL key again.

### Important Note

During calibration, the weighing pan must be unloaded.

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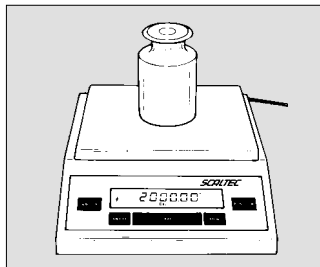
## External Calibration\*\*

Menu code selection: 1 9 1\*

Use only calibration weights with an accuracy equal to or better than the readability of your balance. Refer to the list of calibration weights under "Accessories" on page 73.

When a zero readout is displayed, press the CAL key **(10)**. This starts calibration. The calibration weight value in grams is then displayed.

Errors or interference at the start of the calibration routine are indicated by the error message "Err 02." If this is the case, tare and press the CAL key again when a zero readout appears.

Center the calibration weight on the weighing pan. The balance then calibrates automatically. When calibration is completed, the value of the calibration weight and the stability symbol "g" are displayed.

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\* = factory setting for SBC balances with a readability of  $\leq 0.1$  mg

\*\* = not applicable to verified balances of accuracy class  $\text{II}$

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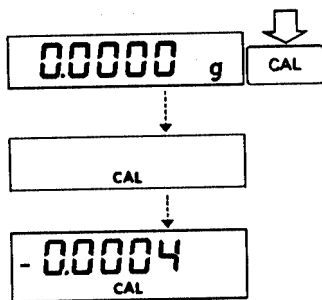
## Calibration Test for SBC Models with a Built-in Calibration Weight

Menu code selection: 1 9 4

Rather substantial changes in barometric pressure and temperature may affect the display response of the balance. To ensure that you obtain the full accuracy of the balance, even when you use the entire weighing range, we have implemented a calibration test function. It will help you quickly to decide whether or not you need to recalibrate your balance (for example, to maintain the same accuracy during long-term weighing series).

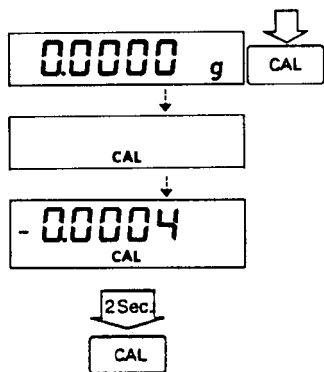
Unload the balance and tare.

When a zero readout is displayed, press the CAL **(10)** key. The built-in calibration weight is now internally applied by servomotor. At the same time, **CAL** is displayed. After the display has stabilized, the deviation of the momentary readout from the target weight (in grams only) is indicated.

If external interference affects the calibration test, you may obtain a brief display of the error message "Err 02." In this case, tare and then press the CAL key again.

### Important Note

The balance should be calibrated if the deviation of the readout from zero is more than the reproducibility specified for your balance in the "Technical Specifications" starting on page 61.

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To calibrate the balance with the built-in weight:  
press the CAL key briefly

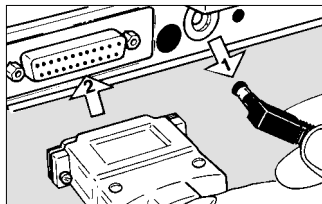
**or**

to quit the calibration test:  
hold down the CAL key for more than 2 seconds

### Blocking the Calibration Functions

You can block the calibration functions by setting code 1 9 7  
(when the menu access switch **(4)** is locked).

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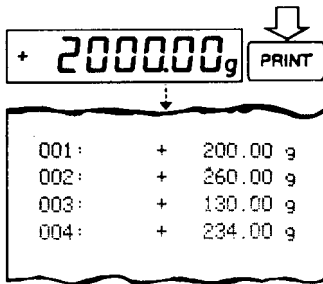
## Data Interface Port

If you wish to record weighing data using a SCALTEC printer, plug the printer connector into the interface port **(15)** of the balance.

### Important Note Concerning Verified Balances Approved for Use as Legal Measuring Instruments:

When using the balance as a legal measuring instrument, you may connect it only to auxiliary measuring devices that are legally permitted. However, if you are not using the balance as a legal measuring instrument, you may connect it to any peripheral devices that you wish, e.g., additional remote displays, personal computer, etc. Such peripheral devices must be marked to indicate that they may not be connected when the balance is being used as a legal measuring instrument.

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### Important Note

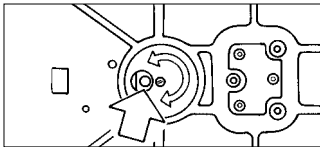
Make sure to unplug the balance from AC power before you connect or disconnect a peripheral device (printer or PC) to or from the interface port.

To print data on hard copy or output them on a computer screen, press the PRINT key **(7)**.

For information about the data output parameters and data ID codes, see page 36. For details on the data interface (such as the data output or input formats, pin assignment, etc.), contact SCALTEC.

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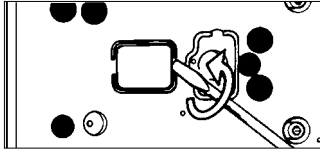
### Below-Balance Weighing

A port for a below-balance weighing hanger is located on the bottom of the balance.

#### For Balances with an Analytical Draft Shield Chamber:


To open the below-balance port, turn the cover plate.

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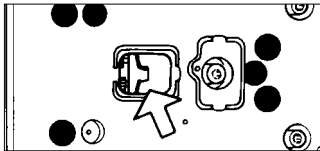**For Balances of the SBA/SBC Series  
(Except for SBA 61/62):**

To open the below-balance port, remove the cover plate from  
the bottom of the balance.

Now you can attach a sample using a suspension wire, for  
example. Common applications for below-balance weighing  
include density determination and immersing a sample  
in a special atmosphere (medium for reaction).

**Important Note Concerning Verified Balances  
Approved for Use as Legal Measuring Instruments:**

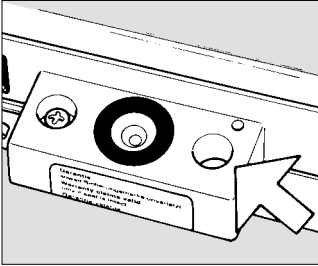
The below-balance weighing port may not be opened or  
used when an approved balance is being operated as a  
legal measuring instrument.

**Important Note**

When you use a below-balance weighing hanger, you must  
install a shield to protect against drafts.

## Diebstahlsicherung\*

## Fastening an Antitheft Locking Device\*

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To fasten an antitheft locking device, use the lug **(17)** on the rear panel of the balance.

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## Troubleshooting Guide

<b>Problem</b>	<b>Causes</b>	<b>Solution</b>
No segments appear on the weight display <b>(11)</b>	<ul style="list-style-type: none"><li>– No AC power is available</li><li>– The AC adapter is not plugged in</li></ul>	<ul style="list-style-type: none"><li>– Check the AC power supply</li><li>– Plug in the AC adapter</li></ul>
No segments appear on the weight display after adjustment/calibration?	<ul style="list-style-type: none"><li>– The surface on which the balance rests is not stable</li><li>– Internal stability has not been reached</li></ul>	<ul style="list-style-type: none"><li>– Make sure that the ambient conditions are stable</li><li>– Prevent vibrations from affecting the surface on which the scale rests</li><li>– Close the draft shield</li></ul>
The weight display shows “H”	<ul style="list-style-type: none"><li>– The load exceeds the capacity of the balance</li></ul>	<ul style="list-style-type: none"><li>– Unload the balance</li></ul>
The weight display shows “L” or “Err 54”	<ul style="list-style-type: none"><li>– The weighing pan <b>(1)</b> and/or the pan support disk is not in place</li></ul>	<ul style="list-style-type: none"><li>– Position the pan and/or the pan support disk (depending on the balance model)</li></ul>
The weight display briefly shows “Err 02”	<ul style="list-style-type: none"><li>– The display did not show a zero readout when the CAL key <b>(10)</b> was pressed to calibrate</li><li>– The balance is loaded</li></ul>	<ul style="list-style-type: none"><li>– Press the TARE key <b>(8)</b>; then press the CAL key again</li><li>– Unload the balance</li></ul>
The special code “◇” does not go out on the weight display	<ul style="list-style-type: none"><li>– None of the keys has been pressed since the balance was turned on</li></ul>	<ul style="list-style-type: none"><li>– Press a key</li></ul>
The weight readout changes constantly	<ul style="list-style-type: none"><li>– Unstable ambient conditions</li><li>– Too much vibration or the balance is exposed to a draft</li></ul>	<ul style="list-style-type: none"><li>– Set up the balance in another area</li><li>– Access the menu to select the correct code for the weighing environment</li></ul>
The weight readout is obviously wrong	<ul style="list-style-type: none"><li>– The balance is not calibrated</li><li>– The balance was not tared before weighing</li><li>– The air bubble of the level indicator <b>(16)</b> is not centered within the circle</li></ul>	<ul style="list-style-type: none"><li>– Calibrate the balance (see page 17)</li><li>– Tare before weighing</li><li>– Level the balance (see page 12)</li></ul>

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### **Servicing**

Regular servicing by a SCALTEC service technician will extend the service life of your balance and increase the measuring accuracy. SCALTEC can offer you service contracts, with your choice of regular maintenance intervals ranging from 1 month to 2 years.

### **Safety Inspection**

If there is any indication that safe operation of the balance with the AC adapter is no longer warranted, turn off the power and disconnect the equipment from AC power immediately. Lock the equipment in a secure place to ensure that it cannot be used for the time being.

In this case, notify your SCALTEC customer service office.

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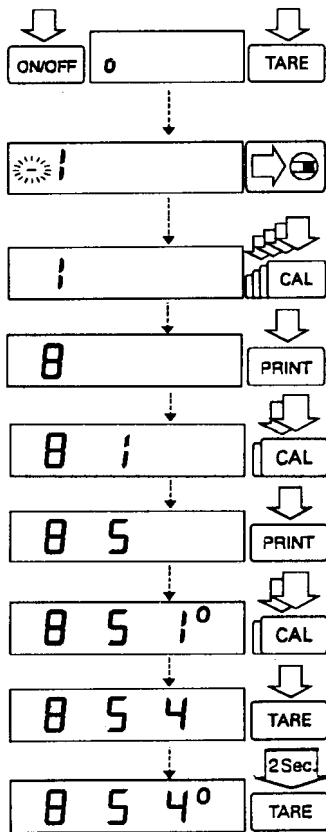
- Accessing the menu
- Setting a code
- Confirming and storing this code

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### Accessing the Menu

- Press ON/OFF to turn off the balance.  
Turn the balance back on. While all segments are displayed, briefly press the TARE key (**8**).
- Release the TARE key when "!" is displayed.
- If "-" is displayed in front of the left-hand number, remove the protective cap from the rear panel of the balance, and move the menu access switch (**4**) in the direction of the arrow.
- Press CAL until "8" appears
- Press PRINT until the 2nd number of the code appears
- Press CAL until "5" appears
- Press PRINT until the 3rd number appears (when you move to this number, the previously set menu code will appear)
- Press CAL to select "4"
- **Confirming a Menu Code Setting**  
Press TARE to confirm the code you have just set (this is indicated by the "o" after the code).
- To **store** the new menu code setting, press TARE for more than 2 seconds.

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### Important Note

The current code setting in the balance operating menu is identified by a small, superscript "°" after the last number. When you access the operating menu, the previously set code will be displayed after you have selected the right-hand number.

If you would like to change several menu code settings, you do not have to press the TARE key after each change to exit the balance operating menu.

Please do not forget to relock the balance operating menu. The symbol "-" indicates that the menu is locked.

### Undoing All Menu Code Changes - Reset Function

The reset function lets you undo all menu code changes, which means that you will obtain the original factory-set menu codes identified by an "\*" To use this function, select code **9 – – 1°**.

Reset Function	Code
On	9 – – 1
Off	9 – – 2*

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<b>Adapting the Balance to Ambient Conditions</b>	Code
Very stable conditions	1 1 1*
Stable conditions	1 1 2*
Unstable conditions	1 1 3
Very unstable conditions	1 1 4

<b>Standard Weighing Mode – Manual Filling Mode</b>	Code
Standard weighing mode	1 2 1*
Manual filling mode	1 2 2

<b>Stability Range</b>	Code
0.25 digit	1 3 1
0.5 digit	1 3 2*
1 digits	1 3 3
2 digits	1 3 4*
4 digits	1 3 5
8 digits <sup>1)</sup>	1 3 6

<b>Tare Parameter</b>	Code**
At any time	1 5 1*/**
Not until the readout is stable	1 5 2*

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\*\* = setting not applicable to verified balances of accuracy class (II)

1) = not applicable to the verified balances approved for use as legal measuring instruments

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<b>Calibration and Linearization Functions Using CAL</b>	Code
External calibration	1 9 1*/**
Internal calibration for SBC models	1 9 3*
Calibration test for SBC models	1 9 4
Internal linearization for SBC 22	1 9 6
Calibration functions blocked	1 9 7

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## Weighing in Two Ranges

The FUNCTION key **(6)** lets you switch back and forth (toggle) between two weight units. The 2nd unit is identified by the display symbol "R1."

	Code
FUNCTION key blocked	2 1 1 *
Mass unit conversion by toggling	2 1 2 *

### Weight Units

The "basic weight unit" is the unit in which your balance will weigh the moment you turn it on. This unit is selected for the first range. You can select a different unit for each of the two ranges by setting the appropriate menu codes.

	Symbol	Conversion Factor 1 g =	1st Range	Code 2nd Range R1
Grams (o)	o	1.	1 7 1	3 1 1
Grams	g	1.	1 7 2 *	3 1 2 *
Kilograms <sup>1)</sup>	kg	0.001	1 7 3	3 1 3
Carats**	ct	5.	1 7 4	3 1 4
Pounds**	lb	0.0022046226	1 7 5	3 1 5
Ounces**	oz	0.035273962	1 7 6	3 1 6
Troy ounces**	ozt	0.032150747	1 7 7	3 1 7
Hong Kong taels**tl		0.02671725	1 7 8	3 1 8
Singapore taels**tl		0.02646063	1 7 9	3 1 9
Taiwanese taels**tl		0.02666666	1 7 10	3 1 10
Grains**	GN	15.43235835	1 7 11	3 1 11
Pennyweights**	dwt	0.643014931	1 7 12	3 1 12

	Symbol	Conversion Factor 1 g =	1st Range	Code 2nd Range R1
Milligrams <sup>2)</sup>	mg	1000.	1 7 13	3 1 13 *
Parts per pound**o		1.1287667712	1 7 14	3 1 14
Chinese taels**tl		0.02645547175	1 7 15	3 1 15
Mommes**	m	0.2667	1 7 16	3 1 16
Austrian carats**k		5.	1 7 17	3 1 17
Tola**	t	0.0857333381	1 7 18	3 1 18
Baht**	b	0.06578947436	1 7 19	3 1 19
Mesghal**	m	0.217	1 7 20	3 1 20

### Display Mode

	Code 1st Range	Code 2nd Range R1
Highest possible accuracy <sup>3)</sup>	1 8 1 *	3 2 1 *
Rounding factor 2**	1 8 3	3 2 3
Rounding factor 5**	1 8 4	3 2 4
Rounding factor 10**	1 8 5	3 2 5

\* = factory setting; depends on the balance model in some cases

\*\* = not applicable to verified balances approved for use as legal measuring instruments

1) = not applicable to verified balances of accuracy class **(I)**

2) = not applicable to verified balances of accuracy class **(II)**

3) = applicable only to verified balances approved for use as legal measuring instruments

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n

<b>Baud Rate</b>	Code
150 baud	5 1 1
300 baud	5 1 2
600 baud	5 1 3
1,200 baud	5 1 4*
2,400 baud	5 1 5
4,800 baud	5 1 6
9,600 baud	5 1 7
19,200 baud	5 1 8
<b>Parity</b>	Code
Mark parity	5 2 1
Space parity	5 2 2
Odd parity	5 2 3*
Even parity	5 2 4
<b>Number of Stop Bits</b>	Code
1 stop bit	5 3 1*
2 stop bits	5 3 2
<b>Handshake Mode</b>	Code
Software handshake	5 4 1
Hardware handshake with 2 characters after CTS	5 4 2*
Hardware handshake with 1 character after CTS	5 4 3

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## Utilities for Printouts or Data Transfer

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<b>Data Output Parameter</b>	Code
Print on request regardless of stability	6 1 1
Print on request after stability, with storage of the function	6 1 2*
Print on request at stability, without storage of the function	6 1 3
Auto print regardless of stability	6 1 4
Auto print at stability	6 1 5
<b>Auto Print</b>	Code
Start/stop auto print using the PRINT key	6 2 1
Auto print not stoppable	6 2 2*
<b>Data Output at Defined Intervals</b>	Code
1 display update	6 3 1*
2 display updates	6 3 2
<b>Automatic Taring after Data Output</b>	Code
Data output without automatic taring	6 4 1*
Data output with automatic taring	6 4 2
<b>Automatic Output of the Application Parameters</b>	Code
Off	7 1 1*
On	7 1 2

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## Additional Functions

### Menu Access Function

You can define the function of the menu access switch by setting the code for the balance operating menu to "accessible." The balance operating menu will then be accessible at all times, which means that you can change the menu codes at any time regardless of the setting of the menu access switch.

<b>Access to Balance Operating Menu</b>	Code
Accessible (codes can be changed)	8 1 1
Locked (codes can be read only)	8 1 2*

<b>Key Functions</b>	Code
Accessible	8 3 1*
Blocked (except for ON/OFF)	8 3 2

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## Menu Access Switch Function on Verified Balances of Accuracy Class Ⅱ

The balance operating menu can also be changed when the balance is being used as a legal measuring instrument. Codes that are not permitted for operation of the balance as a legal measuring instrument are blocked and cannot be selected. The balance operating menu on verified balances approved for use as legal measuring instruments cannot be locked with the menu access switch (“-” not displayed). For balances of accuracy class Ⅱ, the menu access switch is adjusted to lock the “External Calibration” function after verification. To do this, the switch must be moved to the left. A control seal is affixed to the balance to lock the menu access switch. Unauthorized attempts to remove this seal will irreversibly damage it. If you break the seal, the validity of the verification will become void, and you must have your balance re-verified.

Power-on Mode	Code
(Power) off -> on <-> standby	8 5 1 *
on <-> standby	8 5 3
Automatic power-on	8 5 4

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## Application Programs

In addition to the functions implemented for weighing, SCALTEC balances offer you a variety of application programs.

**To select an application program or a specific function, set the appropriate code in the menu.**

Additional settings for running a particular program and displaying or printing the data on hard copy are listed in a table of codes.

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## Tare Memory

### Tare Memory

Press the FUNCTION key **(6)** to store the tare weight. The balance is now automatically tared so you can weigh starting with a zero readout. If you have stored a value in the tare memory, a "1" will be printed or output after the ID code of this value to identify it as such on the printout or on the computer screen.

### "Tare – Net – Gross" Weights

Menu codes used in the example:	Code
Tare memory	2 1 6
Automatic output of all parameters	7 1 2
With data ID code	7 2 2
Print tare memory	7 3 2

As an alternative, you can select this output parameter:

Print net value	7 3 1*
-----------------	--------

### Code 2 1 6

**Application:** Obtaining printouts of tare, net and gross weights

Step/Key	Readout	Data Output/Printout
CF <sup>1)</sup> , TARE	0.00 g	
Place container on pan; press FUNCTION	+ 22.65 g 0.00 g NET	T 1 + 22.65 g
Fill container with sample; press PRINT	+ 250.24 g NET	N 1 + 250.24 g
CF <sup>1)</sup> , PRINT	+ 272.89 g	N + 272.89 g
The data ID codes		mean:
T 1 +	22.65 g	tare weight stored in the memory (weight value)
N 1 +	250.24 g	net weight when tare weight is stored
N +	272.89 g	gross weight = tare + net

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## Practical Example: Net Total

### Menu codes used in the example:

	Code
Tare memory	2 1 6
Automatic data output	7 1 2
With data ID code	7 2 2
Printout of individual components/tare weights	7 3 1*

As an alternative, you can select this output parameter:

Printout of the net total weight/total tare weight	7 3 2
--	-------

**Application:** Simple compounding and formulating of several components with additive storage, automatic taring and outputting of component weight

Step/Key	Readout	Data Output/Printout		
Place container on pan	+ 22.65 g			
CF <sup>1)</sup> , TARE	0.00 g			
Weigh in 1st component; store: FUNCTION	+ 4.61 g			
	0.00 g	NET N 1	+	4.61 g
Weigh in 2nd component; store: FUNCTION	+ 60.33 g			
	0.00 g	NET N 1	+	60.33 g
Weigh in and store additional components: FUNCTION	xxx.xx g	NET N 1	+	xxx.xx g
Finish weighing in components; output total weight: CF <sup>1)</sup> , PRINT	+ 272.89 g	N	+	272.89 g
The data ID codes		mean:		
		N 1	+	60.33 g
		N	+	272.89 g
				component weight (net individual weight)
				total components weighed-in (net total weight)

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## Weighing in Percent

Code 2 1 5

This application program enables you to obtain weight readouts in percent which are in proportion to a reference weight. The reference weight readout is stored as a menu-defined percentage (factory setting: 100%).

### Changing the Reference Percentage

You can change the reference percentage in cycles. Choose from the following settings: 5, 10, 20, 50 and 100.

To activate the change function: Hold down FUNCTION **(6)** for more than 2 seconds until "rEF 100%" appears on the display

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To store the percentage: Hold down FUNCTION for more than 2 seconds to store it permanently (in the non-volatile memory) after you turn off the power

This setting is **not** canceled by the reset code 9 – – 1°!

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With three decimal places

3 6 4

If the weight stored is too light to be displayed, the number of decimal places is automatically decreased.

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3 5 2

**Counting****Code 2 1 4**

The counting program allows automatic conversion of weights into piece counts based on a reference sample weight. A weight readout is stored as a reference sample quantity (factory setting: 10 pcs = pieces).

**Changing the Reference Sample Quantity**

You can change the reference sample quantity in cycles. Choose from the following settings for the reference sample quantity: 5, 10, 20, 50 and 100.

To activate the change function: Hold down FUNCTION **(6)** for more than 2 seconds until "rEF ... pcs" appears on the display

To change the setting: Briefly press FUNCTION again

To store the quantity: Hold down FUNCTION for more than 2 seconds to store it permanently (in the non-volatile memory) after you turn off the power

This setting is **not** canceled by the reset code 9 — 1°!

**Storage Parameter for the Reference Sample Weight**

Code

With full accuracy according to the internal resolution

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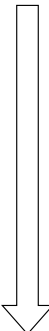
## Practical Example: Parts Counting

### Menu codes used in the example:

	Code
Counting	2 1 4
Reference quantity and ref. weight	7 1 2
With data ID code	7 2 2

**Application:** Counting bulk quantities of items with the same weight

Step/Key	Readout	Data Output/Printout
Place container on pan	+ 22.65 g	
CF <sup>1)</sup> , TARE	0.00 g	
Add 10 counted parts	rEF 10 pcs	

Change the reference sample qty., if desired: To use 20 as the reference qty., add 10 more parts to container. To change the ref. qty. to 20, hold down FUNCTION for >2 seconds rEF 10 pcs

<sup>1)</sup> = press CF for more than two seconds

Press FUNCTION rEF 20 pcs  
To store: hold down FUNCTION for more than 2 seconds

Press FUNCTION to confirm reference piece count + 10 pcs

<b>nRef</b>	+	<b>10 pcs</b>
<b>wRef</b>	+	<b>5.6546 g</b>

Fill container with desired quantity of parts (without counting them) + 500 pcs

<b>PRINT</b>	+	<b>500 pcs</b>	<b>Qnt</b>	+	<b>500 pcs</b>
--------------	---	----------------	------------	---	----------------

Press CF to clear the weight readout and the reference value + 2827.35 g

<b>PRINT</b>	+	<b>2827.35 g</b>
--------------	---	------------------

The data ID codes mean:

Data ID Code	Meaning
<b>nRef</b> + <b>10 pcs</b>	reference piece count
<b>wRef</b> + <b>5.6546 g</b>	reference weight
<b>Qnt</b> + <b>500 pcs</b>	calculated piece count
<b>N</b> + <b>2827.35 g</b>	net weight

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## Animal Weighing/Averaging

Code: 2 1 7

Use this program to determine the weights of live animals or weights under unstable ambient conditions.

In this program, the balance calculates the weight as the average of a selectable number of individual weighing operations by the balance. These weighing operations are also known as "subweighing operations."

During averaging, the number of individual subweighing operations left is shown in the application display in a "countdown" mode.

Once all subweighing operations have been performed, the calculated mean value is indicated as a stable readout on the weight display.

### Manual or Automatic Start Mode

Depending on the menu code you select, animal weighing will be started automatically or manually by pressing a key.

If you select the "automatic mode," you will have to press FUNCTION to start the program for averaging the first weight. You can press CF to interrupt a weight measurement in progress at any time. In the "automatic mode," the symbol "**AUTO**" will be displayed during weight measurement.

The result is locked into the display. The "mouse" symbol or "**AUTO**" flashes during this time. The readout will stop flashing after you have unloaded the balance, and the next weight will be displayed.

### Animal weighing/averaging will start by:

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## Delayed Start Mode

A rule of thumb to go by for selecting the right setting to weigh animals is: the more active an animal is, the greater the difference must be between two successive subweights measured.

Depending on individual requirements, starting the averaging operation can be delayed either in the automatic or manual mode until the animal you are weighing has calmed down to a certain degree.

In this case, the start criterion is defined by the difference between two successive subweights measured.

If the animal moves, the start criterion is not met; therefore, averaging will not start.

Once the animal has calmed down, the program checks whether two measured subweights are within the previously selected range. If so, the actual averaging operation will be started.

Delay start	Code
Slight	3 7 1
Average	3 7 2*
Considerable	3 7 3

During averaging, the number of subweighing operations left to perform is shown on the weight display (countdown mode).

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## Changing the Number of Subweighing Operations

You can change the number of subweighing operations used to average a weight. Change this number in cycles. You can choose from the following settings: 5, 10, 20, 50 and 100.

To activate the change function: Hold down FUNCTION **(6)** for more than 2 seconds until "rEF 10" appears on the display  
If you enter a wrong number, press CF to clear it.

To change the setting: Briefly press FUNCTION

To store the number: Hold down FUNCTION for more than 2 seconds to store this number permanently (in the non-volatile memory) after you turn off the power

This setting is **not** canceled by the reset code 9 — — 1°!

## Storage Threshold for the Automatic Start Mode

To obtain an added measure of reliability in the automatic mode in order to avoid an "incorrect start," a weight value must correspond to a minimum load of 100 display increments.

Once averaging has been completed, the program will stop until the balance is unloaded to half the value (50 display increments) of the storage threshold.

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## Practical Example: Animal Weighing with Automatic Start Mode

### Menu codes used in the example:

	Code
Animal weighing	2 1 7
Delay start until dif. is average	3 7 2*
Automatic start mode	3 8 2*
Automatic output of all parameters	7 1 2
With data ID code	7 2 2

### As an alternative you can select this code:

	Code
Other delay start	3 7 x
Manual start mode	3 8 1

You can change the number of subweighing operations.

If you do not enter a number, averaging will be done based on the number shown on the display (factory setting: 10 subweighing operations).

Once you press FUNCTION the first time, all further weighing operations will be performed automatically after they have met the start criteria.

**Application:** Automatic weighing of animals based on 20 subweighing operations

Step/Key	Readout	Data Output/Printout
Place animal weighing bowl on balance		
CF <sup>1</sup> ), TARE	0.00 g	
Change number of subweighing operations; e.g., set number to 20:		
Hold down FUNCTION for more than 2 seconds	ref 10	
Press FUNCTION	ref 20	
To store: hold down FUNCTION for more than 2 seconds	ref 20	
Place 1st animal in bowl		
FUNCTION	20 465.20 g	m d e f + 20 x - N e t + 4 6 5 . 2 0 g
Remove 1st animal		

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### Application:

Use of the balance in quality assurance systems in compliance with ISO, GLP, GMP, EN, etc.

The balance can record all completed calibration and adjustment operations and print out data in compliance with the requirements of Good Laboratory Practice (GLP). The balance, interfaced with a data printer or a computer, creates a document that records the date, time, serial number and model number, making it possible to clearly trace data to the balance that generated it and the time at which it was generated.

### Select the GLP/GMP-compliant printout or record mode by setting the respective code in the balance operating menu:

GLP/GMP-compliant printout/record mode	Code
Off	8 10 1 *
Only for the calibration and linearization functions	8 10 2
Always on	8 10 3

The following menu code setting must be selected in order to obtain a GLP/GMP-compliant printout/record:

With data ID code	7 2 2
-------------------	-------

### Important Note

GLP/GMP printouts/records will not be generated if the factory setting, code 7 2 1, is selected. In addition, do not select the "Auto print" data output parameter (code 6 1 4 or 6 1 5).

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### **Operating the Balance with a Device that Has a GLP/GMP Printing or Recording Capability:**

You can connect a special SCALTEC printer (order no. SPR 01) to SCALTEC balances for generating GLP/GMP-compliant printouts.

The GLP/GMP-compliant SCALTEC printer provides you with the following functions:

- Turns the GLP/GMP function on or off
- Date/time
- ID no. for identifying a workstation/operator
- Printouts with balance-specific data

### **Printout/Record for Calibration and Linearization Functions**

A printout or record is generated at the end of the following functions:

- All calibration and linearization functions

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The printout can have the following lines:

-----  
**SCALTEC** : Balance family and manufacturer  
**Model SBC22** : Balance model  
**S/N 70500146** : Serial number of the balance  
**Id** : Space for entering a workstation  
or operator ID  
-----

**Date : 21-Jun-97** : Current date  
**Start: 10:05** : Time at which the application started  
**Cal. : Extern** : Calibration mode (in this case,  
"external calibration")  
**Set. : 200.00000 g** : Calibration weight value  
(only for "external calibration")  
**End : 10:05** : End of application  
**Name :** : Space for signature of the operator  
responsible  
-----

Additional information for other calibration modes:

**Cal. : Intern** : Internal calibration  
**Lin. : Intern** : Internal linearization  
**Lin. : Extern** : External linearization  
**Stat.: Complete** : Status comment for calibration  
or linearization functions  
**Cal. : Test** : Calibration test  
**Diff.:- 0.00004 g** : Data measured during  
the calibration test

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## Data Printout/Record

To have a data record printed out, perform the following:

- Press PRINT (after you have turned on the balance or cleared a function by pressing CF) to output the printout heading
- Press PRINT to output additional data
- To end printout generation and recording of data, press CF

Generation of a GLP/GMP-compliant printout or record is also ended when a calibration operation is started.

For application programs, the reference data (parameters) can be included in the printout/record.

<b>Automatic Output of the Reference Data (Parameters)</b>	Code
Off	7 1 1*
Reference %/qty. and reference weight	7 1 2

To have a data record printed out, perform the following:

- Press FUNCTION to output the printout heading and reference data (the reference data will be stored at the same time)

If you input and store new reference values while a GLP/GMP-compliant record is being printed out, the new reference data will be output.

- Press PRINT to output weighing data
- Press CF to end printout generation (this also ends once a calibration or linearization operation has been started)
- Then press CF to clear the reference data for the application programs

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The printout of the data record can have the following lines:

-----  
**SCALTEC** : Balance family and manufacturer  
**Model** **SBC22** : Balance model  
**S/N** **70500146** : Serial number of the balance  
**Id** : Space for entering a workstation  
or operator ID  
-----  
**Date** : **10-Jun-97** : Current date  
**Start** : **11:15** : Time at which the application started  
**Ser.** : : Space for entering  
a project number  
**nRef** + **10 pcs** : Reference data (in this case,  
"counting" – see also page 46)  
**wRef** + **0.13400 g**  
**Qnt** + **500 pcs** : Measured values (in this case,  
"calculated piece count")  
:  
:  
**End** : **12:15** : End of application  
**Name** : : Space for signature of the  
operator responsible  
-----

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## Technical Specifications

Model		SBA 31	SBA 32	SBA 33	SBA 41	SBA 42
Weighing capacity	g	220	150	65	410	120
Readability	g	0.0001	0.0001	0.0001	0.001	0.001
Tare range (subtractive)	g	-220	-150	-65	-410	-120
Repeatability (standard deviation)*	g	≤±0.0001	≤±0.0001	≤±0.0001	≤±0.003	≤±0.003
Linearity	g	≤±0.0002	≤±0.0002	≤±0.0002	≤±0.002	≤±0.002
Response time (average)	s	≤2	≤2	≤2	≤1.5	≤1.5
Adaptation to ambient conditions and application requirements		By selection of 1 of 4 optimized filter levels				
Display update (depends on the filter level selected)	s	0.1 - 0.4				
Allowable ambient operating temperature	°C	+5 ... +40	+5 ... +40	+5 ... +40	0 ... +40	0 ... +40
Operating temperature range	°C	+10 ... +30				
Sensitivity drift within +10 ... +30 °C	/°C	≤±2 · 10 <sup>-6</sup>				
Pan size	mm	Ø 80	Ø 80	Ø 80	Ø 115	Ø 115
Weighing chamber height	mm	225	225	225	-	-
Net weight, approx.	kg	5.4	5.4	5.4	2.8	2.8
Selectable weight units		g, kg, ct, lb, oz, ozt, tlh, tls, tlt, GN, dwt, mg, /lb, tlc, mom, K, tol, bat, MS				
Selectable application programs		Mass unit conversion by toggling; tare memory; net total; weighing in percent; animal weighing				
Automatic zero tracking function (can be turned off by menu code)		Standard feature				
External calibration weight value (of at least accuracy class...)	g	200 (E2)	100 (E2)	50 (E2)	200 (F1)	100 (F1)
Power requirements	V-	Via AC adapter 230 VAC or 115 VAC, -20% ... +15%				
Frequency	Hz	48-63				
Power consumption (average)	VA	16: maximum; 8: average				
Hours of operation with fully charged SBP 01 external battery pack, approx.: h		32				
Built-in interface port		RS-232 C-S/V24-V28; 7-bit; -even, -mark, -odd, -space; Transmission rates: 150 ... 19,200 baud; 1 or 2 stop bits; software/hardware handshake				
<b>Standard features/equipment supplied:</b>						
Analytical draft shield chamber / Dust cover		x	x	x	-	-
Round glass draft shield		-	-	-	x	x
AC adapter, varies acc. to country		x	x	x	x	x
Level indicator		x	x	x	x	x
Hanger for below-balance weighing		x	x	x	x	x

\* = standard deviation of the reproducibility acc. to DIN 8120, Part 3

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## Technical Specifications

Model		SBA 51	SBA 52	SBA 53	SBA 61	SBA 62
Weighing capacity	g	4100	2220	620	7200	4200
Readability	g	0.01	0.01	0.01	0.1	0.1
Tare range (subtractive)	g	-4100	-2220	-620	-7200	-4200
Repeatability (standard deviation)*	g	≤±0.03	≤±0.03	≤±0.03	≤±0.2	≤±0.2
Linearity	g	≤±0.02	≤±0.02	≤±0.01	≤±0.2	≤±0.1
Response time (average)	s	≤1.5	≤1.5	≤1.5	≤1	≤1
Adaptation to ambient conditions and application requirements		By selection of 1 of 4 optimized filter levels				
Display update (depends on the filter level selected)	s	0.1-0.4				
Allowable ambient operating temperature	°C	0...+40				
Operating temperature range	°C	+10...+30				
Sensitivity drift within +10...+30 °C	/°C	≤±2 · 10 <sup>-6</sup>	≤±2 · 10 <sup>-6</sup>	≤±2 · 10 <sup>-6</sup>	≤±5 · 10 <sup>-6</sup>	≤±5 · 10 <sup>-6</sup>
Pan size	mm	180 x 180	180 x 180	Ø 115	180 x 180	180 x 180
Net weight, approx.	kg	3.5	3.5	1.9	2.8	2.8
Selectable weight units		g, kg, ct, lb, oz, ozt, tth, tls, tlt, GN, dwt, mg, /lb, tlc, mom, K, tol, bat, MS				
Selectable application programs		Mass unit conversion by toggling; tare memory; net total; weighing in percent; animal weighing				
Automatic zero tracking function (can be turned off by menu code)		Standard feature				
External calibration weight value (of at least accuracy class...)	g	2000 (F1)	2000 (F1)	500 (F2)	5000 (F2)	2000 (F2)
Power requirements	V~	Via AC adapter 230 VAC or 115 VAC; -20%...+15%				
Frequency	Hz	48-63				
Power consumption (average)	VA	16: maximum; 8: average				
Hours of operation with fully charged SBP 01 external battery pack, approx.:	h	25	25	25	45	45
Built-in interface port		RS-232 C-S/V24-V28; 7-bit; -even, -mark, -odd, -space; Transmission rates: 150...19,200 baud; 1 or 2 stop bits; software/hardware handshake				
<b>Standard features/equipment supplied:</b>						
AC adapter, varies acc. to country		x	x	x	x	x
Level indicator		x	x	x	x	x
Hanger for below-balance weighing		x	x	x	-	-

\* = standard deviation of the reproducibility acc. to DIN 8120, Part 3

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## Technical Specifications

### Verified Balances with EC Type Approval

Model	SBC 22	
Accuracy class		Ⓢ
Weighing capacity Max.*	g	60/210
Scale interval d*	mg	0.01/0.1
Verification scale interval e*	g	0.001
Minimum capacity Min.*	g	0.001
Tare range (subtractive)	g	-210
Range of use according to Directive No. 90/384/EEC*	g	0.001-210
Response time (average)	s	≤12/3
Adaptation to ambient conditions and application requirements	By selection of 1 of 4 optimized filters	
Display update (depends on the filter level selected)	s	0.2-0.4
Allowable ambient operating temperature	°C	+5...+40
Operating temperature range	°C	+15...+25
Pan size	mm	Ø 80
Weighing chamber height (effect. dimens.)	mm	225
Net weight approx.	kg	6,6
Selectable weight units	g, mg	
Selectable application programs	Mass unit conversion by toggling; tare memory; net total; weighing in percent; animal weighing	
Automatic zero tracking function (can be turned off by menu code)	Standard feature	
Power requirements	V-	Via AC adapter 230 VAC or 115 VAC, -20% ... +15%
Frequency	Hz	48-63
Power consumption (average)	VA	16: maximum; 8: average
Hours of operation with fully charged SBP 01 external battery pack, approx.: h	25	
Built-in interface port	RS-232 C-S/V24-V28; 7-bit; -even, -odd; Transmission rates: 150 ... 19200 baud; 1 or 2 stop bits; software/hardware handshake	
<b>Standard features/equipment supplied:</b>		
Dust cover	x	
AC adapter, varies acc. to country	x	
Built-in calibration weight	x	
Level indicator	x	

\* = EC Directive No. 90/384/EEC non-automatic weighing instruments applicable in the E.U. and within the European Economic Area

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## Technical Specifications

Model		SBC 31	SBC 32	SBC 33	SBC 41	SBC 42
Accuracy class		Ⓡ	Ⓡ	Ⓡ	Ⓢ	Ⓢ
Weighing capacity Max.*	g	220	150	65	410	120
Scale interval d*	mg	0.1	0.1	0.1	1	1
Verification scale interval e*	g	0.001	0.001	0.001	0.01	0.01
Minimum capacity Min.*	g	0.01	0.01	0.01	0.02	0.02
Tare range (subtractive)	g	-220	-150	-65	-410	-120
Range of use according to Directive No. 90/384/EEC*	g	0.01-220	0.01-150	0.01-65	0.02-320	0.02-120
Response time (average)	s	≤2	≤2	≤2	≤1.5	≤1.5
Adaptation to ambient conditions and application requirements		By selection of 1 of 4 optimized filters				
Display update (depends on the filter level selected)	s	0.1-0.4				
Allowable ambient operating temperature	°C	+5...+40	+5...+40	+5...+40	0...+40	0...+40
Operating temperature range	°C	+15...+25	+15...+25	+15...+25	+10...+30	+10...+30
Pan size	mm	Ø 80	Ø 80	Ø 80	Ø 115	Ø 115
Weighing chamber height	mm	225	225	225	-	-
Net weight approx.	kg	5.3	5.3	5.3	4.7	4.7
Selectable weight units		g, mg	g, mg	g, mg	g, kg	g, kg
Selectable application programs		Mass unit conversion by toggling; tare memory; net total; weighing in percent; animal weighing				
Automatic zero tracking function (can be turned off by menu code)		Standard feature				
Power requirements	V-	Via AC adapter 230 VAC or 115 VAC, -20% ... +15%				
Frequency	Hz	48-63				
Power consumption (average)	VA	16: maximum; 8: average				
Hours of operation with fully charged SBP 01 external battery pack, approx.: h		32				
Built-in interface port		RS-232 C-S/V24-V28; 7-bit; -even, -odd; Transmission rates: 150 ... 19200 baud; 1 or 2 stop bits; software/hardware handshake				
<b>Standard features/equipment supplied:</b>						
Analytical draft shield		x	x	x	-	-
Dust cover		x	x	x	-	-
Round glass draft shield		-	-	-	x	x
Built-in calibration weight		x	x	x	x	x
AC adapter, varies acc. to country		x	x	x	x	x
Level indicator		x	x	x	x	x

\* = EC Directive No. 90/384/EEC for Non-Automatic Weighing Instruments applicable in the E.U. and within the European Economic Area

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## Technical Specifications

Model	SBC 51	SBC 52	SBC 53	SBC 61
Accuracy class	II	II	II	II
Weighing capacity Max.*	g 4100	2200	620	6200
Scale interval d*	g 0.01	0.01	0.01	0.1
Verification scale interval e*	g 0.1	0.1	0.1	1
Minimum capacity Min.*	g 0.5	0.5	0.5	5
Tare range (subtractive)	g -4100	-2200	-620	-6200
Range of use according to Directive No. 90/384/EEC*	g 0.5-4100	0.5-2200	0.5-620	5-6200
Response time (average)	s ≤1.5	≤1.5	≤1.5	≤1
Adaptation to ambient conditions and application requirements	By selection of 1 of 4 optimized filters			
Display update (depends on the filter level selected)	s 0.1-0.4			
Allowable ambient operating temperature	°C 0...+40			
Operating temperature range	°C +10...+30			
Pan size	mm 180 x 180	180 x 180	Ø 115	180 x 180
Net weight approx.	kg 4.4	4.2	4.7	4.4
Selectable weight units	g, kg			
Selectable application programs	Mass unit conversion by toggling; tare memory; net total; weighing in percent; animal weighing			
Automatic zero tracking function (can be turned off by menu code)	Standard feature			
Power requirements	Via AC adapter 230 VAC or 115 VAC, -20% ... +15%			
Frequency	Hz 48-63			
Power consumption (average)	VA 16: maximum; 8: average			
Hours of operation with fully charged SBP 01 external battery pack, approx.: h	25			
Built-in interface port	RS-232 C-S/V24-V28; 7-bit; -even, -odd; Transmission rates: 150 ... 19200 baud; 1 or 2 stop bits; software/hardware handshake			
<b>Standard features/equipment supplied:</b>				
AC adapter, varies acc. to country	x	x	x	x
Built-in calibration weight	x	x	x	x
Level indicator	x	x	x	x

\* = EC Directive No. 90/384/EEC for Non-Automatic Weighing Instruments applicable in the E.U. and within the European Economic Area

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## Declarations of Conformity

---

**Council Directives of the EU and European Standards:**  
**This SCALTEC balance meets the requirements of the following Directives issued by the Council of the European Union.**

**Council Directive 89/336/EU “Electromagnetic compatibility (EMC)”**

**Important Note:**

The operator shall be responsible for any modifications to SCALTEC equipment and for any connections of cables or equipment not supplied by SCALTEC and must check and, if necessary, correct these modifications and connections. On request, SCALTEC will provide information on the minimum operating specifications (in accordance with the Standards listed above for defined immunity to interference).

**73/23/EEC “Electrical equipment designed for use within certain voltage limits”**

If you use electrical equipment in installations and under ambient conditions requiring higher safety standards, you must comply with the provisions as specified in the applicable regulations for installations in your country.

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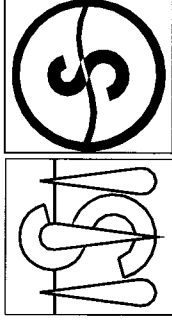
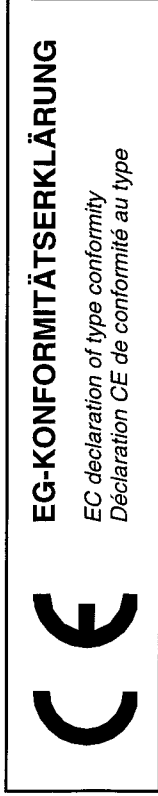
**Additional Directive for Weighing Instruments Used in Legal Metrology: Directive 90/384/EEC “Non-automatic weighing instruments”**

This Directive also regulates the performance of the EC verification by the manufacturer, provided that an EC Type-Approval Certificate has been issued (see page 72) and the manufacturer has been accredited by an officer of a Notified Body registered at the Commission of the European Community for performing such verification. On February 15, 1993, the manufacturers received accreditation for performing verification.

**Subsequent Verifications within the European Countries**

The validity of the verification will become void in accordance with the national regulations of the country in which the balance is used. For information on verification and legal regulations currently applicable in your country, and to obtain the names of the persons to contact, please contact your local SCALTEC office, dealer or service center.

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Copyright by Scaltec Instruments GmbH, Goettingen, Germany.  
All rights reserved.

No part of this publication may be reprinted or translated in  
any form or by any means without the prior written permission  
of Scaltec Instruments GmbH.

The status of the information, specifications and illustrations  
in this manual is indicated by the date given below.

Scaltec Instruments GmbH reserves the right to make changes to the technology,  
features, specifications, and design of the equipment without notice.

Status: March 2001, Scaltec Instruments GmbH, Goettingen, Germany.

**98648-007-63**

Printed in Germany  
on paper that has been bleached  
without any use of chlorine - W198.02 - KT  
Publication No. WXX6032-a01035

**SCALTEC**